



# PocketSwab® Procedural Flow Chart

Refer to Current Charm® Operator's Manual for Complete Test Procedure


Refer to Current Charm® Operator's Manual for Complete Test Procedure

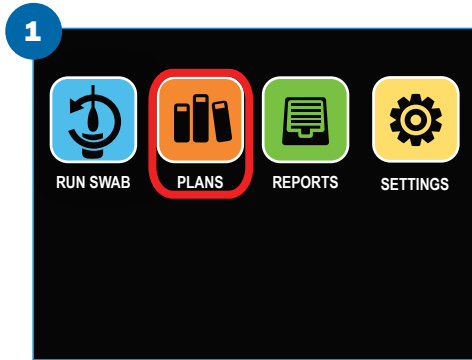


1 Select PLANS button.

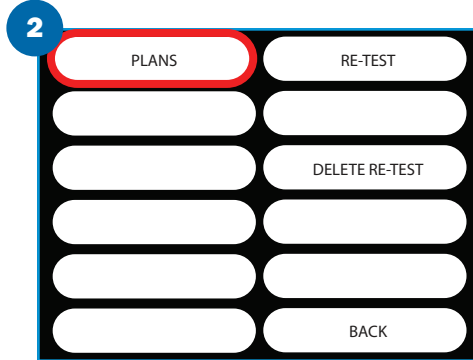

2 Select PLANS.

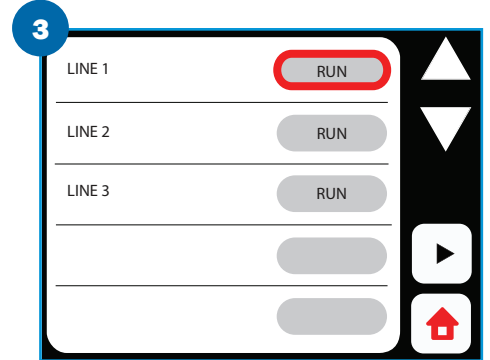

3 Select RUN.

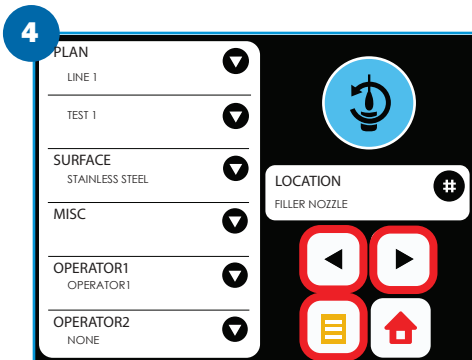

4 Use Left ( ◀ ) or Right ( ▶ ) arrows to pick location or use List View ( ☰ ) to scroll to find correct location.




5 Withdraw the swab by gently pulling and twisting the swab handle.




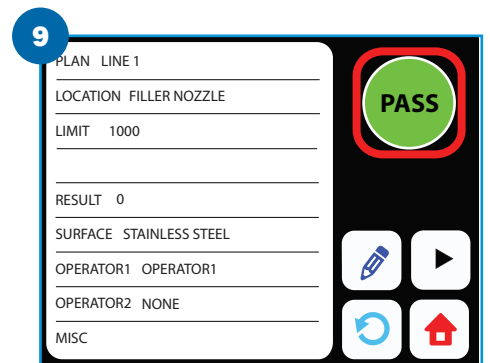
6 Swab an area 100 cm<sup>2</sup> (16 in<sup>2</sup>). Rotate the handle to achieve full swab contact with surface.



7 Reinsert the swab and twist the handle down. Gently shake side to side 3 times. White tablet will break apart.

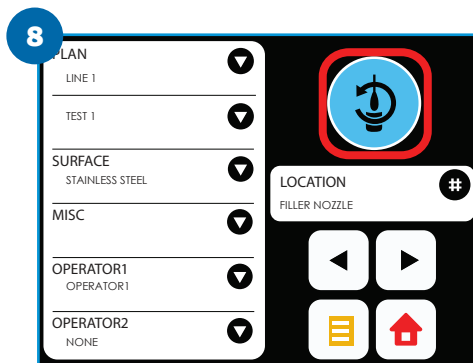

8 Insert the PocketSwab Plus swab fully into the chamber of the novalUM® II system in an upright position. Press  to start count.



9 Results in 5 seconds. The result of "Pass" or "Fail" is based on established limit.

**Note:** Swabs can be held for up to 6 hours at room temperature before activation.



NOTHING WORKS LIKE A CHARM

659 Andover Street, Lawrence, MA 01843-1032, USA  
 T +1.978.687.9200 | F +1.978.687.9216 | E info@charm.com | www.charm.com

© 2016 Charm Sciences, Inc. Charm, novalUM and PocketSwab are registered trademarks of Charm Sciences, Inc.

MRK-1056-001 May-2016